



ELISA for Detection of HTLV-I and HTLV-II Positive Sera

Dr. Michele Owen, HIV Immunology and Diagnostics Branch, Division of AIDS, STD, and TB Laboratory Research, National Center for HIV, STD, and TB Prevention, CDC, 1600 Clifton Rd NE, MS-A25, Atlanta, GA 30333

Reagents

HTLV-I/II Peptides	HTLV-I Peptides Env-1, Env-5, Gag-1a (Catalog# 1459, 1461, 1463); HTLV-II Peptides Env-20, Env-2 (Catalog# 1462, 1460).
Carbonate Buffer	Sigma, Catalog# C3041, dissolve in distilled water.
PBS-T	PBS with 0.05% Tween-20
Blocking Solution	PBS-T with 3% BSA
Antibody Diluent	PBS-T with 1% BSA
Primary Antibody	HTLV-I/II test sera, prepared at 1:20 in antibody diluent.
Secondary Antibody	Goat anti-human IgG(Fc) alkaline phosphatase conjugate (Sigma, Catalog# A9544), prepared at 1:10,000 in antibody diluent.
PNPP Solution	p-Nitrophenyl phosphate (Sigma, Catalog# N2640). Dissolve 1 tablet per 5 ml in sodium carbonate buffer immediately prior to use.
Sodium Carbonate Buffer	1.69 g NaHCO ₃ , 3.51 g Na ₂ CO ₃ , 0.41 g MgCl ₂ (6H ₂ O); add distilled water to bring volume to 2.0 L, adjust pH to 8.6.

Procedure

1. Coat polyvinyl plate (Immulon-II) with HTLV-I or HTLV-II peptides at 10 µg/ml, 50 µl per well, in carbonate buffer. Incubate plate overnight at 4°C.
2. Wash plates six times in PBS-T.
3. Add 150 µl of blocking solution to each well. Incubate plate for 2 hours at room temperature.
4. Wash plates six times in PBS-T.
5. Add 50 µl of diluted serum to each well. Incubate plate overnight at 4°C.
6. Wash plate six times in PBS-T.
7. Add 50 µl of diluted secondary antibody to each well. Incubate plate for 2 hours at room temperature.
8. Wash plate six times with PBS-T.
9. Add 50 µl PNPP to each well and incubate at room temperature.

ALL RECIPIENTS OF THIS MATERIAL MUST COMPLY WITH ALL APPLICABLE BIOLOGICAL, CHEMICAL, AND/OR RADIOCHEMICAL SAFETY STANDARDS INCLUDING SPECIAL PRACTICES, EQUIPMENT, FACILITIES, AND REGULATIONS. NOT FOR USE IN HUMANS.



NIH AIDS Research & Reference Reagent Program
20301 Century Boulevard
Bldg. 6, Suite 200
Germantown, MD 20874
USA

Phone: 240 686-4740
Fax: 301-515-4015
www.aidsreagent.org

-
10. Read the plate at 405 nm as the yellow color approaches an OD of 1.0 in the positive control and test samples. Seropositivity in the test samples is defined as +2 standard deviations over the normal control mean.

ALL RECIPIENTS OF THIS MATERIAL MUST COMPLY WITH ALL APPLICABLE BIOLOGICAL, CHEMICAL, AND/OR RADIOCHEMICAL SAFETY STANDARDS INCLUDING SPECIAL PRACTICES, EQUIPMENT, FACILITIES, AND REGULATIONS. NOT FOR USE IN HUMANS.