

**APPENDIX: CRYOPRESERVATION OF MRA-154**

Note: Only the cryopreservation of immature parasite stage (rings) is viable by this method. We recommend a parasitemia of 3% or higher of ring-stage parasites for cryopreservation.

1. Centrifuge the culture at 1000 × g for 5 minutes.
2. Wash the pellet once with 10 or more volumes of incomplete RPMI 1640 medium. Centrifuge at 1800 × g for 5 minutes and leave sufficient supernatant to resuspend the pellet.
3. To the volume of packed red blood cells, slowly add dropwise one volume of cold (4°C) Glycerolyte 57 solution. Let stand for 5 minutes at room temperature.
4. Add dropwise an additional 4 volumes of cold Glycerolyte 57 solution to the pellet. Mix well and aliquot 0.5 mL into 1.5 mL sterile cryopreservation vials.
5. Place the vials in a controlled-rate freezing unit. From room temperature, cool the vials at -1°C/minute to -40°C. If the freezing unit can compensate for the heat of fusion, maintain rate at -1°C/minute through this phase. At -40°C, plunge vials into liquid nitrogen. Alternatively, place the vials in a Nalgene 1°C freezing container. Place the container at -80°C for 1 to 2 days and then plunge vials into liquid nitrogen.
6. Store in either the vapor or liquid phase of a nitrogen refrigerator (-130°C or colder).